
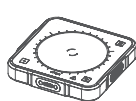
MIXSTER Add-On Unit


1 x MIXSTER
No Power Supply

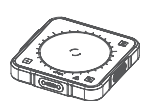MIXSTER Set 1

1 x MIXSTER
1 x Plug-In Power Supply
MIX.PWR.0 (24 W)

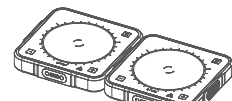MIXSTER Set 2

2 x MIXSTER
1 x Plug-In Power Supply
MIX.PWR.0 (24 W)

MIXSTER Set 3

3 x MIXSTER
1 x Plug-In Power Supply
MIX.PWR.1 (40 W)

MIXSTER Set 4

4 x MIXSTER
1 x Plug-In Power Supply
MIX.PWR.1 (40 W)

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EN Operating instructions Quick Guide

MIXSTER
Attention: This quick guide will inform you about the most important information operations. You can download the complete operating instructions in your language on our website: www.ika.com

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1. EU declaration of conformity

CE We declare under our sole responsibility that this product corresponds to the directives 2014/35/EU, 2006/42/EC, 2014/30/EU and 2011/65/EU and conforms with the following standards or normative documents: EN 61010-1, EN 61010-2-051, EN 61326-1, EN 60529 and EN ISO 12100. A copy of the complete Declaration of Conformity or further declarations of conformity can be requested at sales@ika.com.

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2. Legend of symbols /// Warning Symbols

- DANGER** Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.
- WARNING** Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
- CAUTION** Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
- NOTICE** Indicates practices which, if not avoided, can result in equipment damage.

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3. Assembly

Observe the ambient conditions (temperature, humidity, etc.) listed under "Technical Data".

- Connect the device to a socket.
- Place a magnetic rod in a vessel that is penetrable for magnetic field lines and fill the vessel with the medium to be processed. Then place the vessel on the center of the base plate.

- Notice!**
- Before adding an additional stirring unit, please check your existing power adapter to ensure sufficient power!

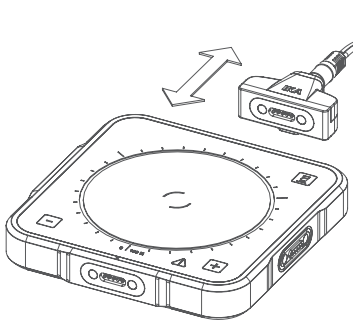
/// Operator panel and display

LED-light: White.....Indicates "stand by" or motor "ON"
LED-light: Green (ON) or Red (OFF).....Indicates "Vessel detection"
LED-light: Blue.....Indicates "setpoint / actual speed"
LED-light: Blue.....Blue-LED blinking = setpoint speed not reached
LED-light: Blue.....Blue-LED lights up constantly = Setpoint speed reached

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
Power ON/OFF

Power Supply connected/disconnected

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
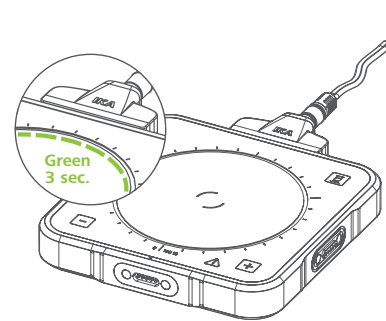
4. MIXSTER Start-up

- Notice!**
- Several device settings can only be changed during the power-up process (LEDs start to light green).
 - The device have no power switch. The power-up process starts as soon as the plug-in power supply unit is connected.

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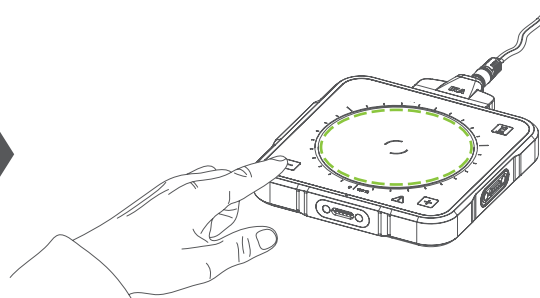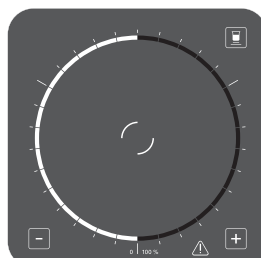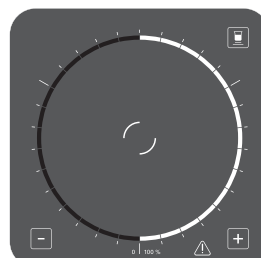
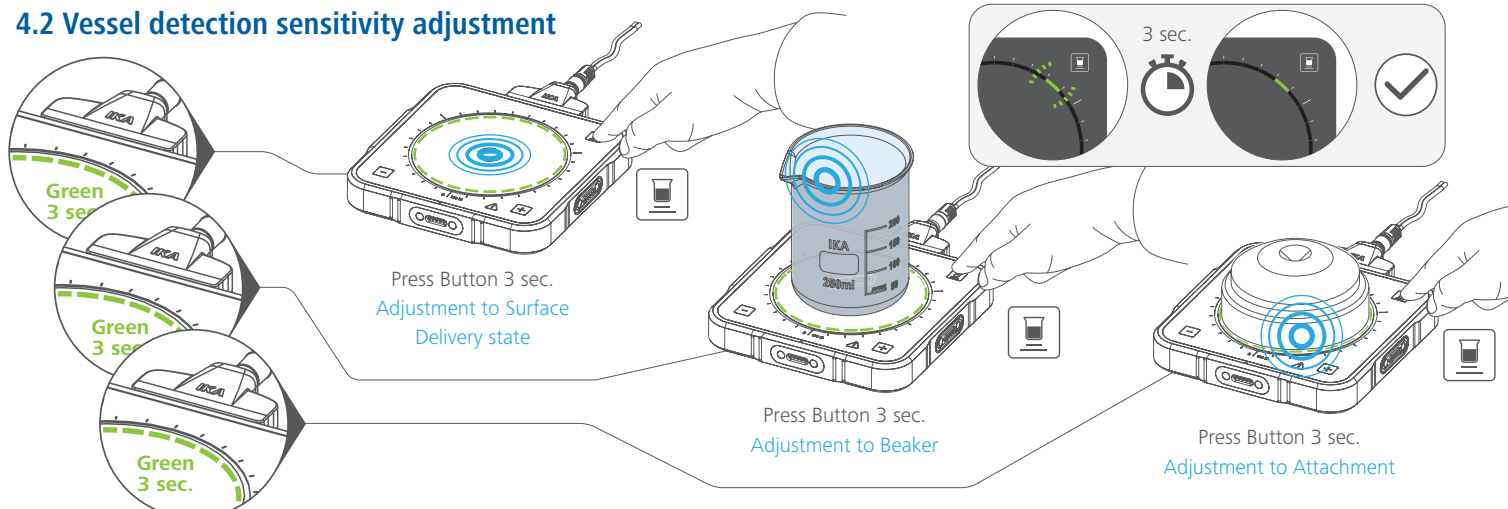
Operating mode A / B

- Mode A
Settings are not saved when the device gets disconnected from the power supply. Stirring function will be set to OFF when the device is powered on.
- Mode B
All settings will be stored if the device gets disconnected from the power supply. The stirring function will be set to ON (if the set value > 0 rpm) or OFF (if the set value = 0 rpm) when the device is powered on, depending on the previous status of the device.

4.2 Vessel detection sensitivity adjustment

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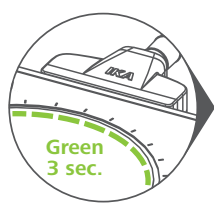
Vessel detection sensitivity adjustment


General setting routine to be performed at the moment of switch-on.
Press and hold the "Vessel detection" button until the green LED stops flashing.

- 4.2.1 Reset or Adjustment to working plate:**
Remove vessel or attachment from the working plate and make the adjustment.
- 4.2.2 Adjustment to a Vessel:**
Place the vessel on the working plate and make the adjustment.
- 4.2.3 Adjustment to an attachment:**
Place the attachment on the working plate and make the adjustment.

4.3 Automatic reverse rotation AR

1 AR ON/OFF Setting

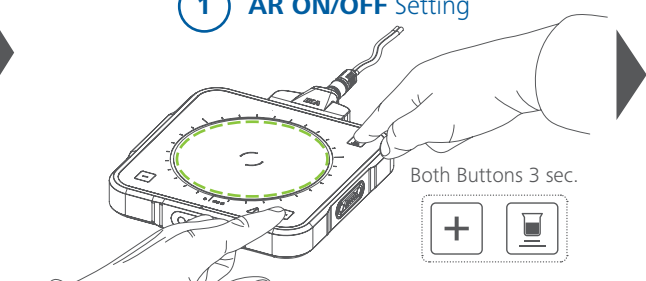Reverse Adjustment Activate

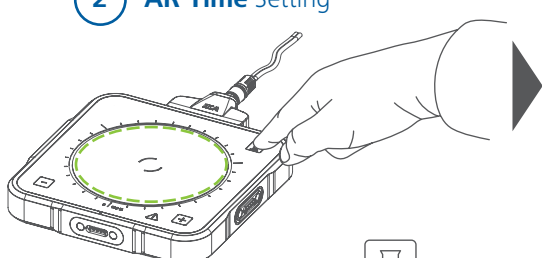
Reverse OFF

max 3 Sec.


Reverse ON

2 AR Time Setting

Timer Activate

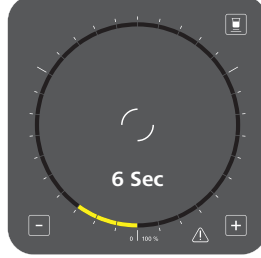Time Setting
Yellow 2 Sec each Digit
Green 1 Min each Digit

Setting Sample
2 Min + 18 Sec
Max Adjustment 30 Min

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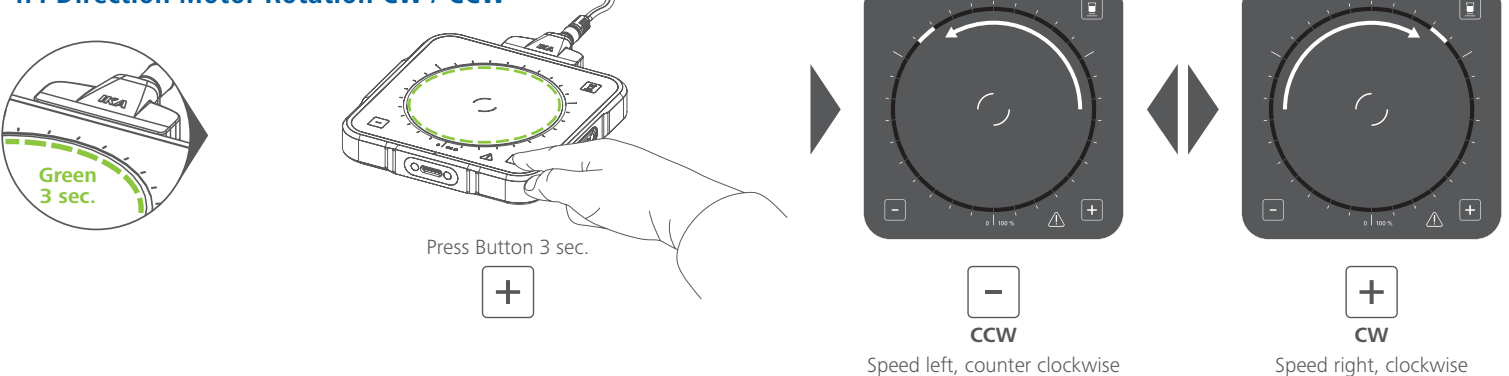
Automatic Reverse AR

Reverse rotation ON / OFF with interval time setting
Continue setting the interval time, after the reverse rotation function has been activated.

4.4 Direction Motor Rotation CW / CCW

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
Rechtsdrehend (CW) / linksdrehend (CCW)



Direction of the Motor Rotation

Clockwise (CW) / counterclockwise (CCW)

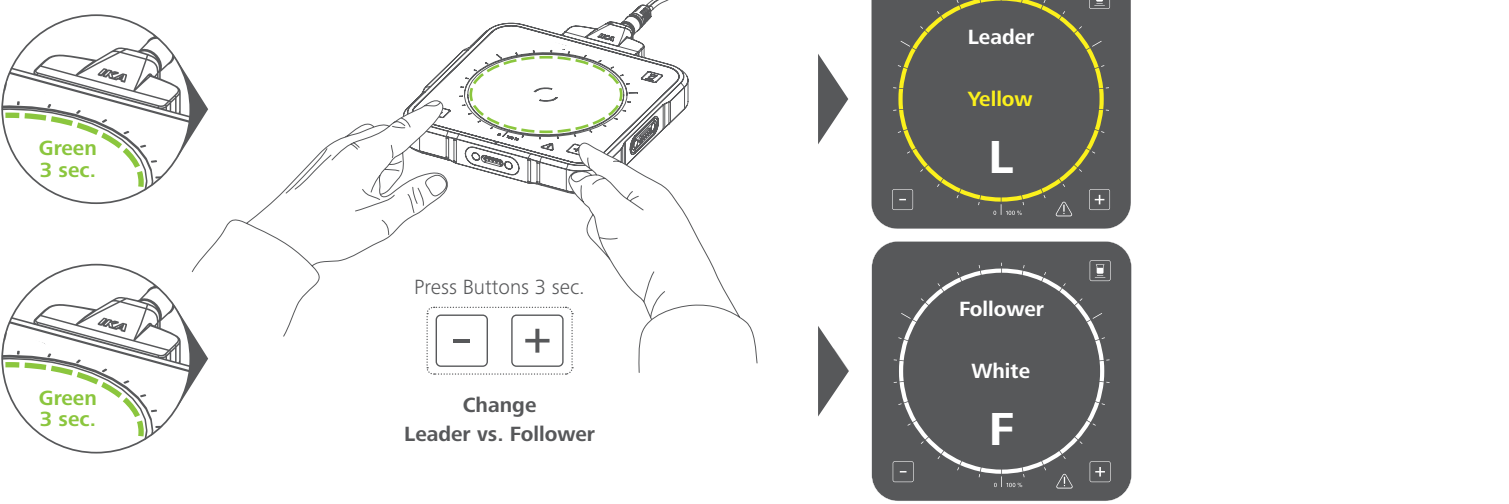4.5 Multi-Position Setup

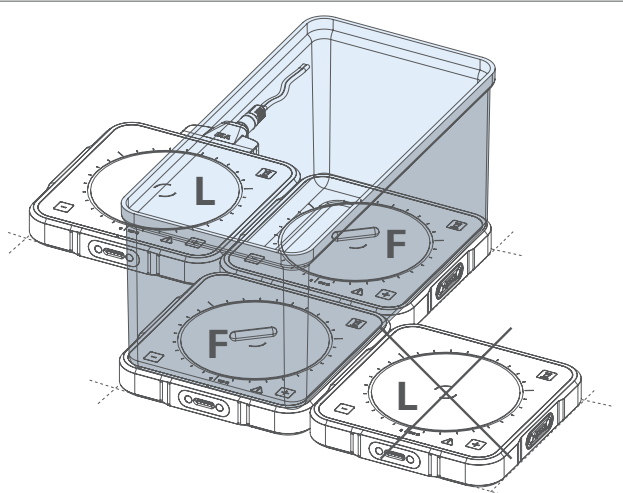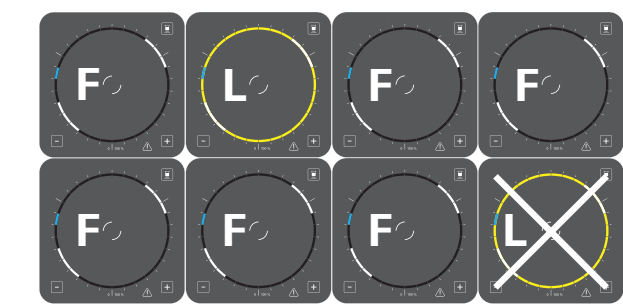NOTICE: Only 1 Leader each multiple arrangement

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Device operation „Follower“ / „Leader“

/// Operation in „Leader“ / „Follower“ mode

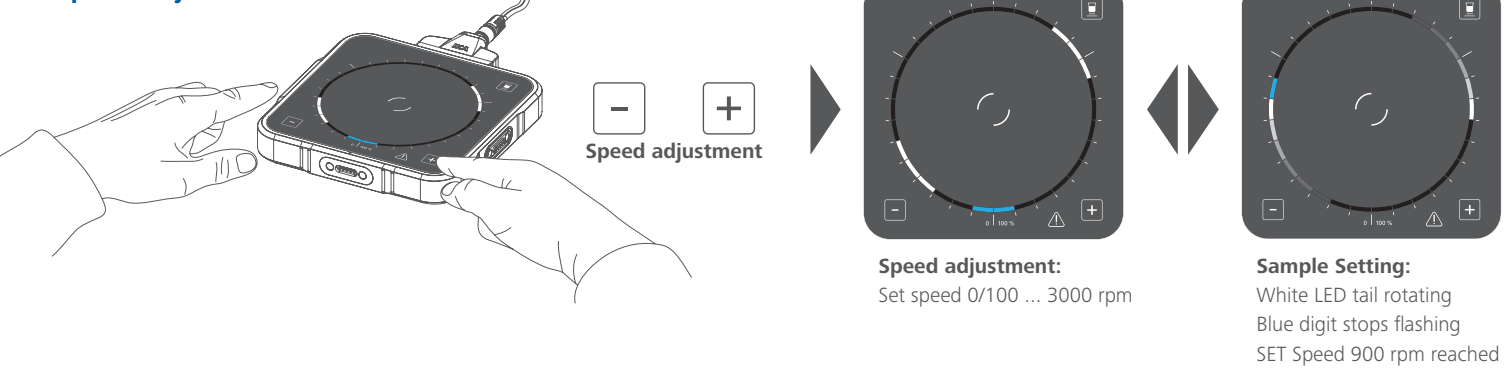
- Notice!**
 - If more than 1 „Leader“ device is detected at the moment of switching-on, all the devices configured as „Leaders“ will be reset to factory settings.
 - Operation of multiple „Leader“ units in one system can damage the device!
 - „Follower“ devices take over the settings of the „Leader“ units.
- Notice!**
 - Once a „Leader“ device has been configured in the stirring system, the „Follower“ devices are controlled synchronously. All keys including the vessel detection of the „Follower“ devices are deactivated.
 - This allows stirring operation of the „Follower“ devices across the permissible footprint, e.g. in the form of a larger container covering the working surface of the „Follower“ devices.

/// Operation in „Leader“ / „Follower“ mode: (L + F) > 4 x

- Notice!**
 - If more than 4 devices [(L + F) > 4 x] are operated at the same time, a more powerful desktop power supply and an additional connector (adapter) are required. (see „Accessories“ www.ika.com.)

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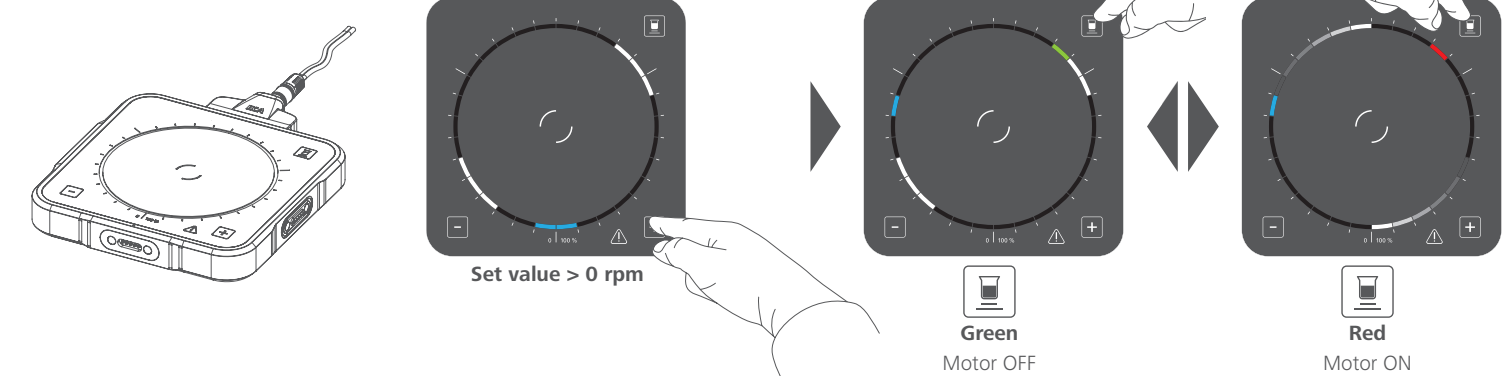
Stirring


- Motor ON: White LEDs with tail, rotate
- Motor OFF: White LEDs, do not rotate (no digit tail)

5.2 Activation of stirring function Following scenarios shows different ways to start the motor function


5.2.1 Activation by key button

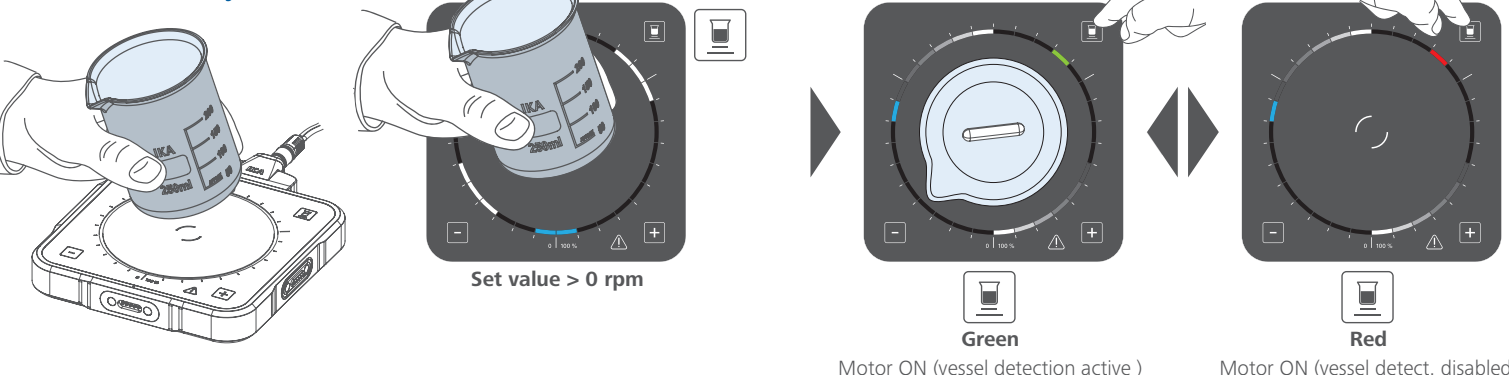
5.2.2 Activation by „vessel detection“

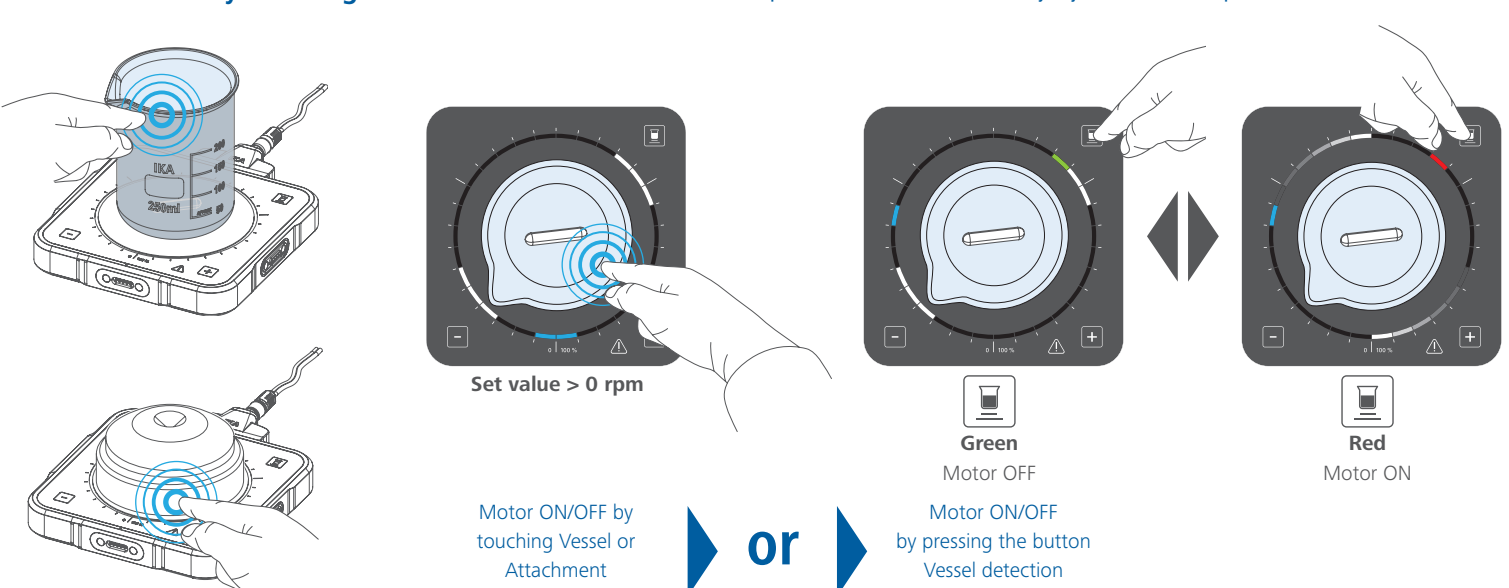5.2.3 Activation by TOUCHING the vessel or the attachment Requires „vessel detection“ sensitivity adjustment > see chapters 4.2.2 and 4.2.3

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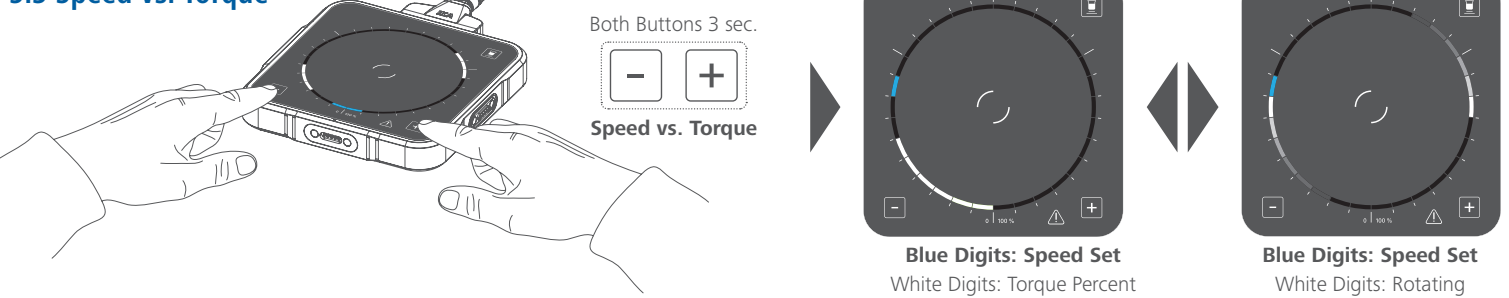Speed vs. Torque

- Selection of speed or torque display
- Relative representation torque. Increase/decrease torques
- Relative representation stirring function. LEDs rotate.

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Maintenance and cleaning

- The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.
- Cleaning: See www.ika.com
- Spare parts order: See www.ika.com
- Repair: See www.ika.com

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Warning / Errors

- Motor overload
 - Maximum internal device temperature exceeded
 - Permitted minimum input voltage is too low
 - Maximum input current allowed overshoot
 - Motor error
- › See www.ika.com

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Technical data

› For detailed data description see www.ika.com

Nominal Voltage	24V DC
Max Power IN	10 W
Motor rating output	5 W
Set-up plate material	Glass
Set-up plate dimension	124 x 124 mm
Dimensions	135 x 135 x 24 mm
Weight	600 g
Permissible ambient temperature	+5 ... +40 °C
Protection Class DIN EN 60525	IP 66

Stirring quantity max H ₂ O	5 L
Speed range	0/100 ... 3000 rpm
Stir bar length	10 ... 30 mm

Power supply input	100 ... 240 VAC, 50/60 Hz
Power supply output	24 VDC, 40/24 W

IKA-Werke GmbH & Co. KG
Janke & Kunkel-Str. 10,
79219 Staufen, Germany
Phone: +49 7633 831-0
eMail: sales@ika.de

USA
IKA Works, Inc.
Phone: +1 910 452-7059
eMail: sales@ika.net

KOREA
IKA Korea Ltd.
Phone: +82 2 2136 6800
eMail: sales-lab@ika.kr

BRAZIL
IKA Brasil
Phone: +55 19 3772 9600
eMail: sales@ika.net.br

VIETNAM
IKA Vietnam Co. Ltd.
Phone: +84 28 38202142
eMail: info_japan@ika.net

JAPAN
IKA Japan K.K.
Phone: +81 6 6730 6781
eMail: info_japan@ika.net.jp

TURKEY
IKA Turkey A.Ş.
Phone: +90 216 394 43 43
eMail: sales.turkey@ika.com

KENYA
IKA Works Kenya Ltd.
Phone: +254 112 323 745
eMail: sales.kenya@ika.com

SPAIN
IKA Works Spain, S. L. Barcelona
Phone: +34 93 6099-5666
eMail: sales.lab@ika.com

MALAYSIA
IKA Works (Asia) Sdn Bhd
Phone: +60 3 6099-5666
eMail: sales.lab@ika.my

CHINA
IKA Works Guangzhou
Phone: +86 20 8222 6771
eMail: info@ika.cn

POLAND
IKA Poland Sp. z o.o.
Phone: +48 22 201 99 79
eMail: sales.poland@ika.com

THAILAND
IKA Works (Thailand) Co. Ltd.
Phone: +66 2059 4690
eMail: sales-lab-thailand@ika.com

INDIA
IKA India Private Limited
Phone: +91 80 26253 900
eMail: info@ika.in

UNITED KINGDOM
IKA England LTD.
Phone: +44 1865 986 162
eMail: sales.England@ika.com

UGANDA
IKA Works Kampala Limited
Phone: +254 112 323 745
eMail: sales.uganda@ika.com

Technical specifications may be changed without prior notice. Discover and order the fascinating products of IKA online: www.ika.com