# Safe Mixing

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\_\_\_\_\_  
 Unterschrift, Firmenstempel

# Index

- About this document ..... 22
  - Typographic conventions ..... 22
  - Copyright protection..... 22
- Basic product information ..... 23
  - Standards and directives, product certification ..... 23
  - Residual risk..... 23
  - Intended use..... 23
  - Reasonably foreseeable misuse..... 23
- Transportation..... 24
- Storage..... 24
- Acclimatization ..... 24
- Permissible ambient conditions ..... 24
- General safety information..... 25
- Electrical safety ..... 25
- Operational safety ..... 25
- Work safety ..... 26
- Personal protective equipment (PPE) ..... 26
- Environmental protection..... 26
- Biohazard ..... 27
- Special hygiene measures for the use of laboratory equipment in food,  
cosmetics and pharmaceutical production ..... 27
  - General Measures ..... 27
  - Device-specific Measures ..... 28
- California Residents ..... 28
- Other regulations ..... 28
- Device overview ..... 29
- Commissioning..... 30
  - Set up the device..... 30
  - Power Supply ..... 30
  - Switch the device on/off..... 30
- Operation ..... 31
- Troubleshooting..... 33

Service work.....	33
Cleaning instructions .....	33
Repairs .....	34
Maintenance.....	34
Disposal .....	34
Technical Specifications .....	35
Scope of delivery .....	35
Accessories.....	35
Contact information Heidolph international.....	36
Warranty Statement .....	36
Declaration of no objection .....	37
Declaration of Conformity .....	38
China RoHS Certification.....	39

## About this document

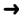
This operating manual describes the features and operation of magnetic stirrers type Hei-MIX S.

The operating instructions manual is an integral part of the delivery!

If applicable, variants will be explicitly identified and described in detail.

## Typographic conventions

Standardized symbols and signal words are used in this document to identify warnings, cautions, and important information.

Symbol	Signal word / explanatory note
	Meanings of the safety symbols and signal words used: <b>DANGER</b> Indicates a hazardous situation which, if not avoided, will result in death or serious injury.
	<b>WARNING</b> Indicates a possible hazardous situation which, if not avoided, could result in death or serious injury.
	<b>CAUTION</b> Indicates a hazardous situation which, if not avoided, could result in minor or moderate injury or material/environmental damage.
	Prohibition signs are used to prevent or avoid a certain activity or situation which could result in injury or material/environmental damage.
	Mandatory signs are used to indicate important information regarding the product handling, This information is used to ensure operational safety and to maintain the value of the product.
	The arrow symbol indicates instructions to be followed in order to ensure the operational safety when handling the product.

## Copyright protection

This publication is protected by copyright and intended for internal use by the purchaser of the product only.

No part of this publication may be transmitted or reproduced in any form, by any means, without the prior written consent of the copyright owner Heidolph Instruments GmbH & Co. KG. Any violation is subject to compensation for damages.

## Basic product information

### Standards and directives, product certification

#### CE Marking

The device complies with the following standards:

- European Machinery Directive, 2006/42/EC
- EMC Directive, 2004/30/EC

### Residual risk

The device was designed and manufactured in accordance with the latest technical standards at the time of development and the recognized safety regulations. During installation and use, as well as during maintenance work, repairs and cleaning, there are nevertheless certain residual risks associated with the device described.

These are identified and described at the appropriate points in this document.

### Intended use

The device described in this manual has been developed and designed by the manufacturer for the following laboratory activities:

- Mixing
- Stirring
- Titrating

Due to its design, the device in its delivery condition may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use of this device is not considered as intended!

### Reasonably foreseeable misuse

For use under conditions or for purposes deviating from the intended use, additional measures may become necessary, and/or specific guidelines and safety regulations will have to be observed (see section "Special hygiene measures for the use of laboratory equipment in food, cosmetics and pharmaceutical production" on page 27). Corresponding requirements must be evaluated and observed by the operator in each individual case.

Compliance with and implementation of all relevant guidelines and safety measures for the respective field of application is within the sole responsibility of the operator.

All risks resulting from improper use are solely borne by the operator.

The device may exclusively be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is operated with responsibility are the sole responsibility of the operator!

## Transportation

During transport, avoid severe shocks and mechanical stresses that can cause damage to the device.

Keep the original packaging in a dry and protected place for later use.

## Storage

Always store the device in its original packaging. To protect against damage and unreasonable material aging, store the device in a dry environment that should be as temperature-stable and dust-free as possible.

Recommended ambient conditions for storage:

- 5 °C – 31 °C up to 80 % rel. humidity
- 32 °C – 40 °C up to 50 % rel. humidity, decreasing linearly

## Acclimatization

After each transport and after storage under critical climatic conditions (high temperature difference between inside and outside), allow the device to acclimatize at room temperature for about two hours to prevent possible damage from condensation before putting it into operation at the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections (power supply, tubing) only after the device has been acclimatized!

## Permissible ambient conditions

The device is designed for indoor use only. Permissible ambient conditions for operation:

- 5 °C – 31 °C up to 80 % rel. humidity
- 32 °C – 40 °C up to 50 % rel. humidity, decreasing linearly
- Maximum height above sea level: 2,000 m

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

The device **IS NOT** suitable for outdoor use!  
The device **IS NOT** suitable for use in hazardous areas!

## General safety information

Observe the following general safety instructions to prevent any personal injury and/or damage to property when handling the device.

- Before commissioning and using the device, familiarize yourself with all the safety regulations and guidelines for occupational safety applicable at the place of use and observe them at all times.
- Only operate the device if it is in perfect technical condition. In particular, make sure that no visible damage can be detected.
- If there is missing or misleading information on the device or on occupational safety, contact the responsible safety specialist or our technical service.
- Only use the device in accordance with the regulations for intended use ("Intended use" on page 23).

## Electrical safety

- Ensure that the voltage indicated on the rating plate matches the supply voltage of the country in which the device is being used.
- Ensure that the mains socket-outlet is protected by means of a residual-current device (RCD).
- Always use the supplied power supply cord provided with the device.
- Prior to use, check that the device and the power supply cord are free of visible damage.
- Have repairs and/or maintenance work on the device carried out exclusively by an authorized and skilled electrician or by the technical service department of Heidolph Instruments.
- Always switch the device OFF and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

## Operational safety

- Do not make any unauthorized changes or modifications to the device!
- Only use genuine spare parts and accessories, or those expressly approved by the manufacturer!
- Observe all safety instructions and hazard warnings as well as all (handling) instructions contained in this document!

Noncompliance will invalidate any warranty against Heidolph Instruments.

The operator is solely liable for all damage resulting from unauthorized changes or modifications to the unit, from the use of unauthorized or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

## Work safety

- Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.
- Persons with pacemakers must maintain a minimum distance of 10 cm from the device to prevent possible interference.
- Depending on the effort and duration, holding the sample in automatic mode can lead to overloads in the hand and arm area. When using the device, observe the specifications of the EC directive 2002-44 on the permissible vibration load.
- Do not operate any other devices in the immediate vicinity of the device,
  - which can generate electromagnetic fields in the frequency range between  $9 \times 10^3$  Hz and  $3 \times 10^{11}$  Hz,
  - which generate emission or radiation sources in the frequency range  $3 \times 10^{11}$  Hz to  $3 \times 10^{15}$  Hz (in the optical spectral range wavelengths from 1,000  $\mu\text{m}$  to 0.1  $\mu\text{m}$ ),
  - which generate ultrasonic or ionizing waves.
- Do not process substances that could release energy in an uncontrolled manner (e.g. self-ignition).
- Do not process substances in which the energy input through mixing poses a danger.
- Do not work with an open flame in the immediate vicinity of the device (risk of explosion).
- Wipe off any fluid that may have spilled on the device immediately.
- Always switch off the device when it is not in use.

Rectify disturbances or faults on the device immediately.  
 Shut down the device and disconnect it from the power supply if it is not possible to eliminate the disturbance or rectify the fault.  
 In such a case, contact the responsible technical service department.

## Personal protective equipment (PPE)

The operator must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is solely within the operator's responsibility.

## Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid risks to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the responsible personnel is the sole responsibility of the operator!

## Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking of the device with the biohazard warning symbol:

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the responsible personnel is the sole responsibility of the operator!

## Special hygiene measures for the use of laboratory equipment in food, cosmetics and pharmaceutical production

When laboratory equipment is used in the production processes of the food, cosmetics or pharmaceutical industry, special hygiene measures must be taken by the user to avoid sample contamination and to minimize any risk to humans and the environment as far as possible.

Please observe the following recommendations:

### General Measures

- Ensure a clean working and storage environment when handling substances and materials.
- Train all employees in the field of occupational hygiene, document all training measures and check the implementation of all required hygiene measures during operation regularly.
- Use a hygiene control concept such as HACCP (Hazard Analysis and critical Control points). The HACCP comprises the following criteria:
  - Hazard analysis
  - Identification of critical control points
  - Definition of critical limit values
  - Establishment of a system for monitoring and controlling critical hazard control points (CCP)
  - Corrective actions for uncontrollable CCP
  - Establishment of a system to verify the implementation of all HACCP measures
  - Establishment of a system for documenting all associated procedures and protocols

The evaluation of the applicability of the mentioned rules and regulations is within the sole responsibility of the operator!

## Device-specific Measures

- Make sure that even products that are intended for single use only are of sufficient purity.
- Avoid contamination by handling contaminated vessels, apparatus or aids with care.

### Contact information

For further information, please contact our after sales service at any time.

Phone: +49-9122-9920-0

Mail: [sales@heidolph.de](mailto:sales@heidolph.de)

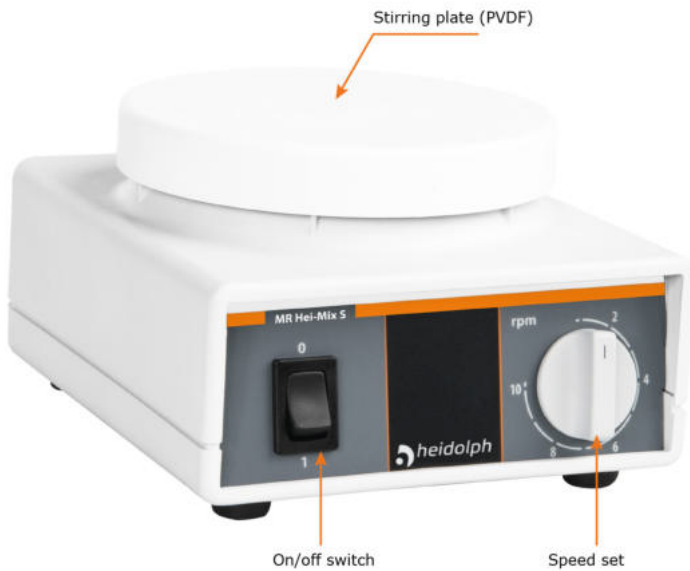
## California Residents

Important information for California residents regarding Prop 65. Please visit [www. P65Warnings.ca.gov](http://www.P65Warnings.ca.gov) for more information.

## Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

## Device overview

## Commissioning

### Set up the device

#### **CAUTION**

#### **Risk Of Slipping!**

The vibration during operation can cause the device to move over the mounting surface and/or fall to the ground.

- Observe the instructions for the correct positioning of the device!

- Place the device on a solid level and horizontal surface for use.
- Before switching on, make sure that the device is sufficiently stable.
- Keep all contact surfaces clean and dry.

### Power Supply

Always use the supplied power supply cord provided with the device! Observe the instructions in section "Electrical safety" on page 25.

The device must only be supplied with power from a properly grounded mains socket-outlet.

### Switch the device on/off

- Set the speed to zero before switching on the device. To do this, turn the speed controller fully anti-clockwise.
- Use the on/off switch to turn the device on or off.

## Operation

Observe the instructions in this section when operating the device!

### CAUTION

#### **Risk of injury, risk of damage to property due to tilting/falling of the device**

High shaking frequencies in combination with a large load and/or a tall attachment assembly on the shaking platform may cause the device to swing up and fall!

- Only gradually increase the shaking frequency to the required level under the above conditions, paying attention to the stability of the structure.
- Lower the shaking frequency or reduce the total load on the shaking platform when the device starts to move during operation.

An uneven distribution of the load on the shaking platform may cause the device to swing up and fall!

- Always ensure that the vessels are evenly distributed on the shaking platform.
- Pay particular attention to an even distribution of the weight load when vessels of different sizes and/or differently filled vessels are placed on the shaking platform at the same time!

#### **Risk of injury, risk of damage to property from falling vessels**

At high shaking frequencies, there is a risk that improperly fastened vessels will fall off the shaking platform.

- Before switching on the device, make sure that all vessels are properly fixed on the shaking platform.

#### **Hand-arm over-exposure to vibration**

Depending on the effort and the duration of use, there is a risk of over-exposure to vibration of your hand and arm when using the hand to hold the vessel during operation.

- Observe the European Directive 2002/44/EC regarding the permissible hand-arm vibration exposure.

**WARNING**

**Risk of injury, risk of damage to property caused by splashing fluids**

When using open vessels, there is a risk that fluid will spray out at high shaking frequencies.

→ Whenever possible, use closed vessels for processing corrosive, toxic or biohazardous substances and seal them safely.

→ Increase the speed only gradually to the desired shaking frequency, especially with open vessels, and observe the fluid movements.

→ Always use the necessary and appropriate personal protective equipment!

**Automatic restart**

When the mains is restored after a power interruption, the device automatically starts up again with the set parameters!

Depending on the setting and structure, there is a risk that the device will move uncontrolled and/or fluid will be sprayed out of the vessels.

→ In the event of an interruption to the mains supply, switch off the device via the main switch.

- Always place individual vessels centrally on the stirrer plate.
- Use suitable attachments for simultaneous processing of several samples/vessels (see section "Accessories" on page 35).
- Do not switch on the unit until all vessels including stirring bars are securely positioned on the stirrer plate.
- Use the speed selector to set the required intensity.

Due to the phase angle control of the unit, the stirring bars in the vessel do not move until the speed selector is at least in position 4.

## Troubleshooting

The following table includes possible failures and corresponding corrective measures.

Failure	Cause	Possible remedy
No stirring function	→ Power supply failure	→ Check power supply cord for correct connection → Check fuse
	→ No stirring bars in the vessel(s)	→ Insert stirring bars
	→ Speed set to 0	→ Increase speed

If a fault cannot be rectified with the described remedies, please contact an authorized sales representative or our technical service (see section "Contact information Heidolph international" on page 36).

## Service work

When carrying out service work on the device (cleaning, maintenance, repair), observe the general instructions and safety information described in this section.

### **WARNING: Danger of electric shock**

Live components are installed inside the device.

When opening the device, there is a risk of touching live components.

- Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

Penetrating liquid poses the danger of an electric shock.

- When cleaning, avoid the penetration of liquids.

## Cleaning instructions

Wipe all surfaces of the device with a damp cloth if necessary. Persistent contamination can be removed with mild soapy water.

### **CAUTION: Damage to the device**

Improper cleaning can damage the surfaces of the device.

Penetrating liquid can damage the electronic components inside the device.

- Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.
- Never use any aggressive or abrasive cleaning agents or aids.

## Repairs

Repairs to the device may only be carried out by authorized skilled experts!

Unauthorized repairs during the warranty period will result in the loss of the warranty claim.

The owner is solely liable for damage caused by unauthorized repairs.

In case of repair contact an authorized dealer or our technical service, see "Contact information Heidolph international" on page 36.

Include the completed declaration of no objection with every device return, see "Declaration of no objection" on page 37.

## Maintenance

The device contains no user-serviceable components. The installed motors are maintenance-free. If necessary, in the event of abnormal operating behavior such as excessive noise or heat generation, for example, contact our technical service.

## Disposal

- When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.
- When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their transposition into national law in the country of use.
- Check the device and all components for residues of substances that are hazardous to health, the environment, and biohazardous before disposing.
- Properly remove and dispose residues of substances that are hazardous to health, the environment and biohazardous!

## Technical Specifications

Supply voltage variants	230 V 50 Hz
	115 V 60 Hz
Power input	7 W
Protection class (IEC 61140)	2 
Protection class (IEC 60529)	IP 21
Acoustic pressure (dB(A))	< 70 (in accordance with IEC 61010)
Drive	shaded pole motor
Operating mode	continuous
Speed range	0 – 2.200 rpm
Speed accuracy	±5 %
Max. stirring capacity (H <sub>2</sub> O)	5 l
Max. load	6 kg
∅ stirrer plate	104 mm
Weight	1.1 kg
Dimensions (W × H × D)	126 × 80 × 140 mm

## Scope of delivery

Component / Variants	Quantity	Product number
<b>Hei-Mix S</b>	1	503-02000-00
Operating instructions English / German	1	01-005-004-60
Guarantee registration / Declaration of no objection	1	01-006-002-78
EC Declaration of Conformity	1	01-001-025-06
Power supply cord	1	country specific

\* The product numbers shown are for 230 V EU devices. For product numbers of variants, please contact Heidolph instruments.

## Accessories

Accessories can be found on our homepage at [www.heidolph-instruments.com](http://www.heidolph-instruments.com)

## Contact information Heidolph international

### Heidolph Instruments North America

Phone: 1-866-650-9604  
E-mail: [service@heidolph.com](mailto:service@heidolph.com)  
[www.heidolphNA.com](http://www.heidolphNA.com)

### Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20  
E-mail: [service@radleys.co.uk](mailto:service@radleys.co.uk)  
[www.heidolph-instruments.co.uk](http://www.heidolph-instruments.co.uk)

### Local distributors

To find your local distributor please visit [www.heidolph.com](http://www.heidolph.com)

## Warranty Statement

Heidolph Instruments assumes a three-year warranty against material and manufacturing defects.

Excluded from the warranty are glass and wear parts, transport damage, and damage resulting from improper handling or non-intended use of the product.

For registered products, the warranty period begins on the date of purchase. Register the product with the enclosed warranty card or on our homepage [www.heidolph.com](http://www.heidolph.com).

For non-registered products, the warranty period begins with the date of the serial production (to be determined by the serial number).

In the event of material or manufacturing defects, the product will either be repaired or replaced free of charge within the warranty period.

## Declaration of no objection

Enclose the declaration of no objection, duly completed, with your device return.  
Submissions without a declaration of no objection cannot be processed!

### 1. Information on the device

Item number: \_\_\_\_\_  
 Serial number: \_\_\_\_\_  
 Reason for submission: \_\_\_\_\_

### 2. Does this product pose a risk to people and/or the environment due to its use for processing substances that are hazardous to health, the environment and/or biohazardous? Please mark with a cross and complete the information!

		If <b>YES</b> , with which substances did the device come into contact?
		_____
		_____
<b>NO</b>	<b>YES</b>	
		If <b>NO</b> , what cleaning and/or decontamination measures were carried out?
		_____
		_____

### 3. Information on the client/sender:

Name, first name: \_\_\_\_\_  
 Company/institution: \_\_\_\_\_  
 Department/working group: \_\_\_\_\_  
 Address: \_\_\_\_\_  
 Zip code, city: \_\_\_\_\_  
 Country: \_\_\_\_\_  
 Phone: \_\_\_\_\_  
 E-Mail: \_\_\_\_\_

### 4. Legally binding declaration

With his signature, the client/submitter declares the completeness and correctness of his details. Missing or incorrect information obligates to compensation for damages.

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hereby declare, that the product (from serial number on 200066828) designated below is in compliance with the basic requirements of all applicable EU-directives stated below with regard to design, type of model sold and manufactured by us. This certificate will be invalid if the product is modified without the prior written consent and agreement of the manufacturer.

MR Hei-Mix S	503-02000-00, ...-01, ...-13, ...-14
MR Hei-Mix L	505-00000-00, ...-13
MR Hei-Mix D	505-01000-00, ...-03
MR Hei-Standard	505-20000-00, ...-02, ...-03, ...-09
MR Hei-Tec	505-30000-00, ...-03
MR Hei-Connect	505-40000-00, ...-16
MR Hei-End	505-50000-00, ...-02, ...-03, ...-05

Carousel Standard RR91200, .../SWISS, .../EURO  
Carousel Tech RR91203, .../SWISS, .../EURO  
Carousel Connect Stirring Hotplate CAT No:RR91207

Maschinenrichtlinie / Machinery Directive 2006/42/EG  
EMV-Richtlinie / Electromagnetic Compatibility Directive 2014/30/EU  
Delegierte (RoHS-) Richtlinie / Delegated (RoHS) Directive 2015/863/EU  
Angewandte (harmonisierte) Normen / (Harmonized) Standards applied:  
EN ISO 12100:2010, EN 61326-1:2013, EN 61010-1:2010, EN IEC 63000:2018

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## China RoHS Certification

### China RoHS DECLARATION OF CONFORMITY

Heidolph Instruments GmbH & Co.KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

#### Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for Heidolph Instruments GmbH & Co.KG products is 25 years.

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This table is created according to SJ/T 11364-2014

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**Note:** Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
- O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
- X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (CrVI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by Heidolph Instruments GmbH & Co.KG may enter into further devices or can be used together with other appliances.

With these products and appliances in particular, Heidolph Instruments GmbH & Co.KG will not take responsibility for the EFUP of those products and appliances.

Place, date Schwabach, 19.09. 2019

Wolfgang Jaenicke  
Chief Executive Officer CEO

Marcell Sarré  
Vice President  
Quality Management &  
Technical Service

© Heidolph Instruments GmbH & Co. KG  
Doc-ID: 01-005-004-60-2 – Ed.: 2021-Aug.-05

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Subject to change without notice. The printed version of this document is not regularly updated. The latest issue of this document can be found by visiting our homepage.

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MATERIAL CONTENT DECLARATION FOR Heidolph Instruments GmbH & Co. KG PRODUCTS